


Compound Library Handling Instructions

Centrifuge

- Products must be thawed and centrifuged before use, because the compound will stick to the tube wall or lid due to bumps and other reasons during transportation. Centrifugation helps reduce unnecessary waste and loss.

Sample pickup

- Our sample rack comes with a laser etched barcode and LidLock on one side. The LidLock can prevent the tubes in racks from falling out if the rack is inverted or dropped. Push the LidLock in the desired direction as marked by the arrows to open or lock the rack.

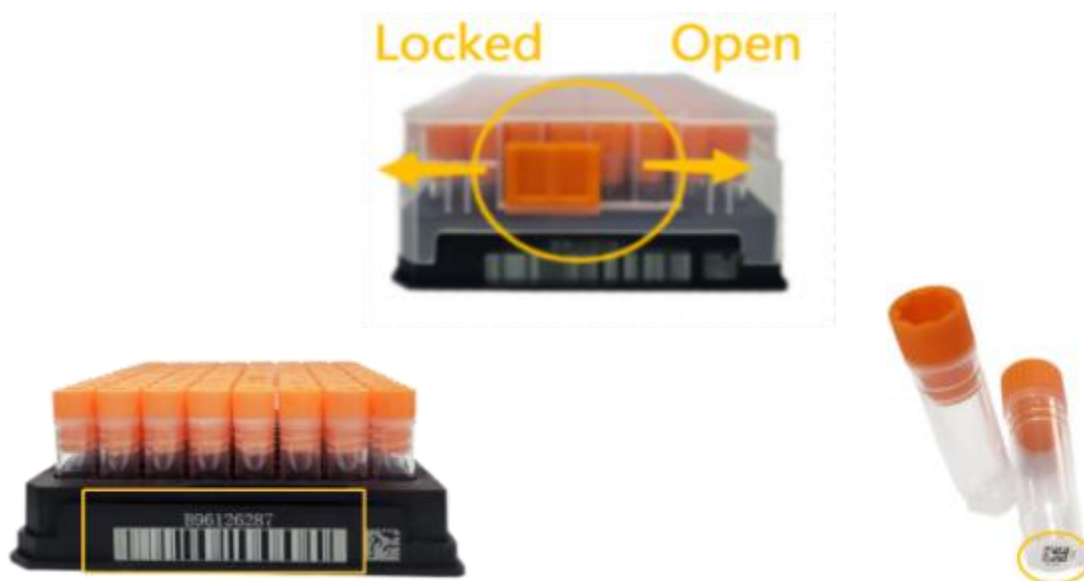

- At the bottom of the freezer tube, there is a unique unrepetitive digital code and the corresponding QR code.

Storage

- Products must be sealed and stored at -20°C or -80°C (Recommended -80°C)
- According to your experimental needs, take part of the mother liquor and dilute it into a sample for experimental operation. The remaining mother liquor and samples shall be sealed and stored at -20°C or -80°C