

Eco-Friendly Wax Remover and Degreaser

(Spray Bottle)

E1507835

Storage: Room Temperature.

Introduction:

The Eco-Friendly Wax Remover and Degreaser can be used to clean residual paraffin and dirt from the surfaces of microtomes, embedding machines, workbenches, and other equipment. Its efficacy is superior to that of imported products. This product does not contain toxic or harmful reagents such as xylene. It can effectively remove paraffin and dirt, and also protect microtomes to extend the service life of the equipment. This reagent is for research use only and not intended for clinical diagnosis or other purposes.

Usage Method (for reference only):

1. Use a pre-cooled medical scraper to scrape off large pieces of paraffin. Use a dental probe to assist in cleaning the gaps, and then use a brush to sweep away the paraffin shavings.
2. Spray this product directly on the surfaces of microtomes, embedding machines, workbenches, and other areas with residual paraffin and dirt. Wait for 5-10 minutes, then wipe clean with toilet paper.

Precautions:

1. This product has a certain degree of volatility. Please use the reagent as soon as possible after opening.
2. For your safety and health, please wear a lab coat and disposable gloves during operation.