

o-Nitrobenzyl Alcohol-modified Sodium Hyaluronate (HANB)

Cat. No.: N1518262 | Pack size: 1 g, 5 g | Storage: Protected from light; store at -20°C

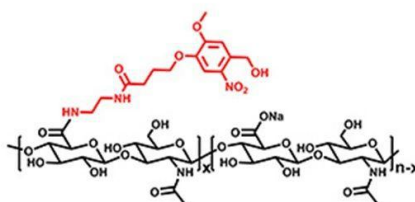


Fig. 1 Chemical structure of o-Nitrobenzyl Alcohol-modified Sodium Hyaluronate (HANB).

Product Introduction

Hyaluronic acid (HA) is an important component of the human extracellular matrix, with key physiological functions including water retention, maintenance of extracellular space, osmotic pressure regulation, lubrication, and promotion of cell repair. HA has been approved by the FDA, exhibits excellent biosafety, and is widely used in the biomedical field.

HANB is a photosensitive polymer material obtained by modifying sodium hyaluronate with o-nitrobenzyl alcohol (NB). Under light irradiation at 365–405 nm, NB undergoes photolysis to produce aldehyde groups that can react with amino groups, thereby enabling covalent conjugation of HANB with amino-containing substances (e.g., proteins, peptides, drugs, bioactive factors) as well as seamless integration and adhesion to tissues. In addition, when used in combination with photosensitive polymers containing double bonds, the photolysis products of NB can participate in free-radical polymerization to form hydrogels with excellent mechanical strength.

Product Specification

Component	Appearance	Specification	Remarks
-----------	------------	---------------	---------

Component	Appearance	Specification	Remarks
HANB	Pale yellow to yellow powder or granules	1 g/vial, 5 g/vial	Store in the dark.

Recommended Use

- HANB can be used in combination with amino-containing polymers such as polylysine, chitosan, carboxymethyl chitosan, and gelatin, and crosslinks to form a gel under light irradiation at 365-405 nm.
- HANB can be used with methacrylate-modified polymers (including CMCMA, GelMA, HAMA, CSMA, ChMA, CMChMA) and allyl-modified polymers such as F127DA, and polymerized under photoinitiator and light exposure to construct hydrogel materials with diverse physicochemical properties.

Preparation Method

- When used with double-bond-containing polymers and photoinitiators, the recommended photoinitiator dosage is:
Low substitution degree HANB (1.5%-2.0%): 30 mg (LAP/NAP) per 1 g HANB;
Medium substitution degree HANB (3.0%-4.0%): 54 mg (LAP/NAP) per 1 g HANB.
- Preparation of photoinitiator stock solution (if applicable): Dissolve an appropriate amount of NAP or LAP to the desired concentration; store at 2-8 °C protected from light.
- Weigh the required amount of HANB into a centrifuge tube, then add the corresponding volume of initiator stock solution or mixture. If clumps form, crush gently with a spatula first.
- Stir the mixture at room temperature in the dark or shake on an orbital shaker until completely dissolved.

Precautions

- Higher substitution degree of HANB and higher solution concentration lead to stronger tissue adhesion, but also higher light blocking and longer required irradiation time.
HANB with substitution degree 1.5%-2.5%: recommended concentration ≤ 10 wt%;
HANB with substitution degree 3.0%-4.0%: recommended concentration ≤ 5 wt%.
- Higher molecular weight of HANB results in higher solution viscosity.
HANB with molecular weight 890 kDa: recommended concentration ≤ 3 wt%;
HANB with molecular weight 340 kDa: recommended concentration ≤ 5 wt%.

Applications

3D chondrocyte culture, hemostatic materials, bone defect repair, drug loading, tissue engineering, regenerative medicine, protein immobilization, bioprinting, postoperative anti-adhesion, etc.

Sterilization Methods

1. Filter sterilization (recommended): Sterilize the solution using a 0.22 μm sterile syringe filter. Only suitable for low-viscosity solutions. For high-viscosity solutions, use pasteurization or autoclaving.
2. Pasteurization: Heat the solution to 80 $^{\circ}\text{C}$ for 30 min, then rapidly cool to room temperature in an ice-water bath. Repeat the cycle three times.
3. Moist heat sterilization (autoclaving): Sterilize in an autoclave at 121 $^{\circ}\text{C}$ for 8 min. After manual venting, cool rapidly to room temperature in an ice-water bath. Note that high temperature may reduce gel performance; strictly follow the above procedure.

Note: Sterile solutions: store at 2-8 $^{\circ}\text{C}$ protected from light, use within 7 days; Non-sterile solutions: store at 2-8 $^{\circ}\text{C}$ protected from light, use within 48 hours.

Product Features

1. In situ formation of high-strength hydrogels: The aldehyde groups generated by HANB photolysis react with amino groups on polymer backbones to form a gel network.

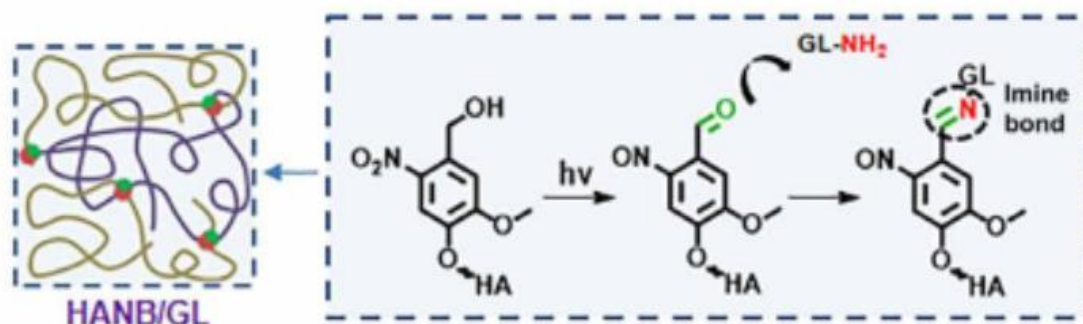


Fig. 2 In situ formation of high-strength hydrogels.

2. Tissue adhesion and integration: Aldehyde groups from HANB photolysis react with amino groups on tissue surfaces at the application site, achieving covalent bonding between the gel and tissue.

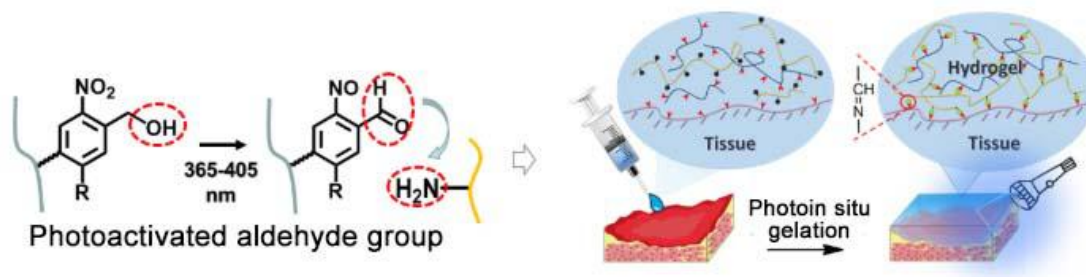


Fig. 3 Tissue adhesion and integration.

3. Immobilization of bioactive substances: Aldehyde groups from HANB photolysis conjugate with amino groups on bioactive molecules, enabling controllable immobilization within the hydrogel.


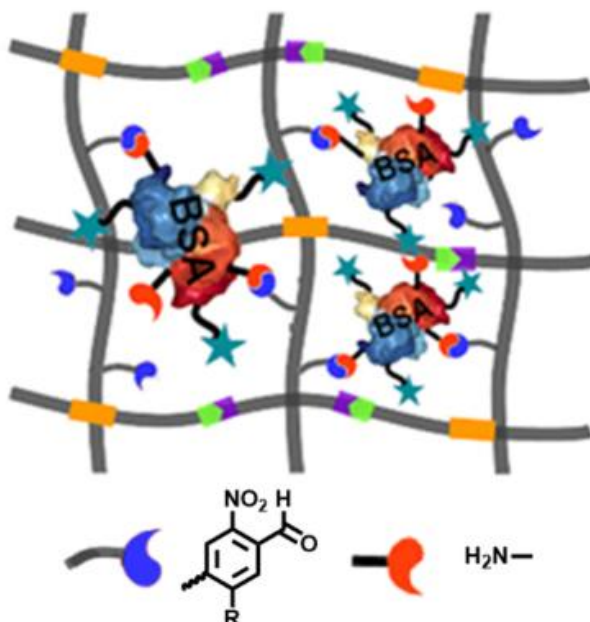


Fig. 4 Immobilization of bioactive substances.

Typical Applications

Benefiting from the above advantages, HANB hydrogel systems have been used by researchers worldwide in cardiac hemostasis, bone/cartilage repair, wound treatment, 3D cell culture, etc., with numerous landmark achievements.


1. Puncture hemostasis: Second-scale curing, high-strength, high-adhesion photocrosslinkable hydrogel for hemostasis of cardiac punctures.

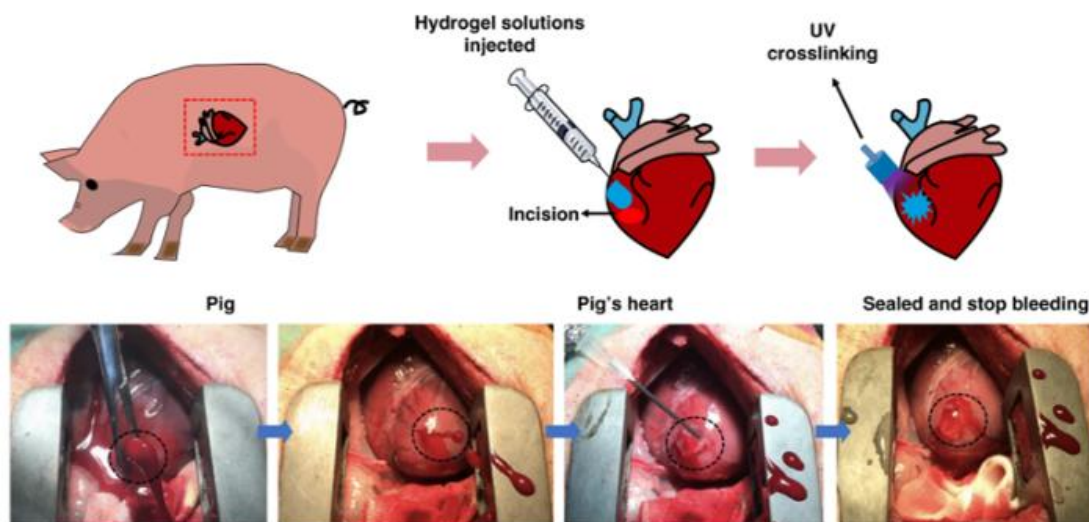

Nat. Commun. (IF = 16.6), 2019, 10, 2060.

Fig. 5 Puncture hemostasis application.


2. Bone repair: Rapid preparation of high-strength cell-laden microgels via microfluidics for bone defect repair.

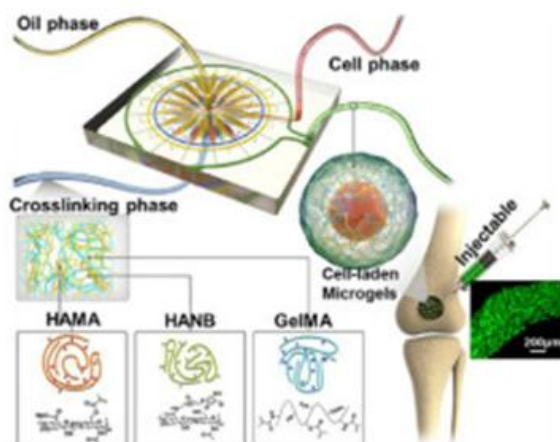

Acta biomaterialia (IF = 9.7), 2023, 157, p.91-107

Fig. 6 Bone repair application.


3. Cartilage repair: Second-scale curing, high-strength, high-adhesion photocrosslinkable hydrogel for arthroscopic cartilage repair.

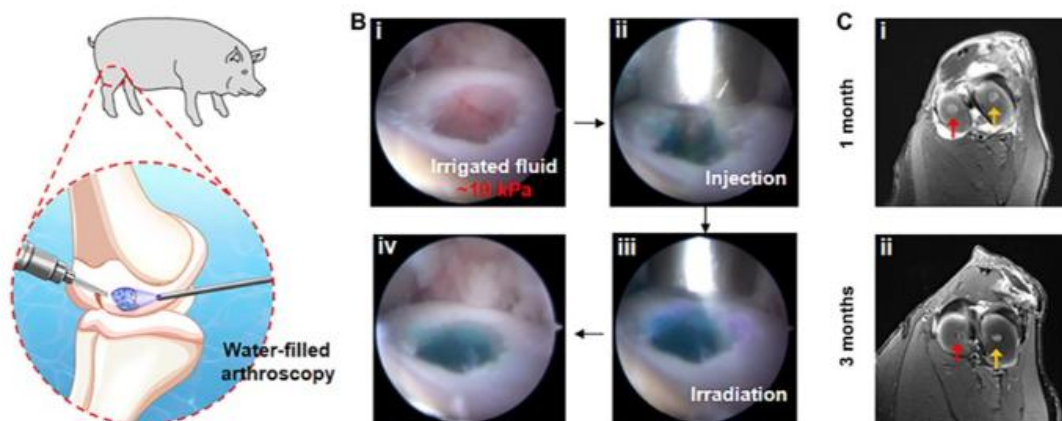

***Sci. Adv.* (IF = 13.6), 2021, 7, eabg0628;**

Fig. 7 Cartilage repair application.

4. Wound treatment: Fast-curing, high-adhesion photocrosslinkable hydrogel for defective wound repair.

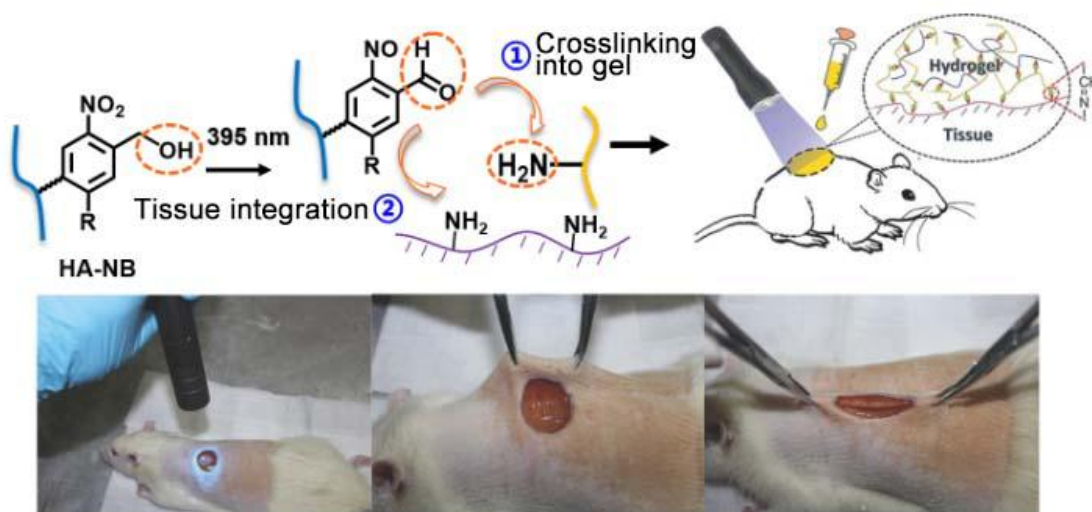

***Adv. Mater.*(IF = 29.4), 2016, 28, 2724.**

Fig. 8 Wound treatment application.

5. Cell culture: Regulation of bioactive molecule immobilization in 3D hydrogels to construct a favorable cellular microenvironment.

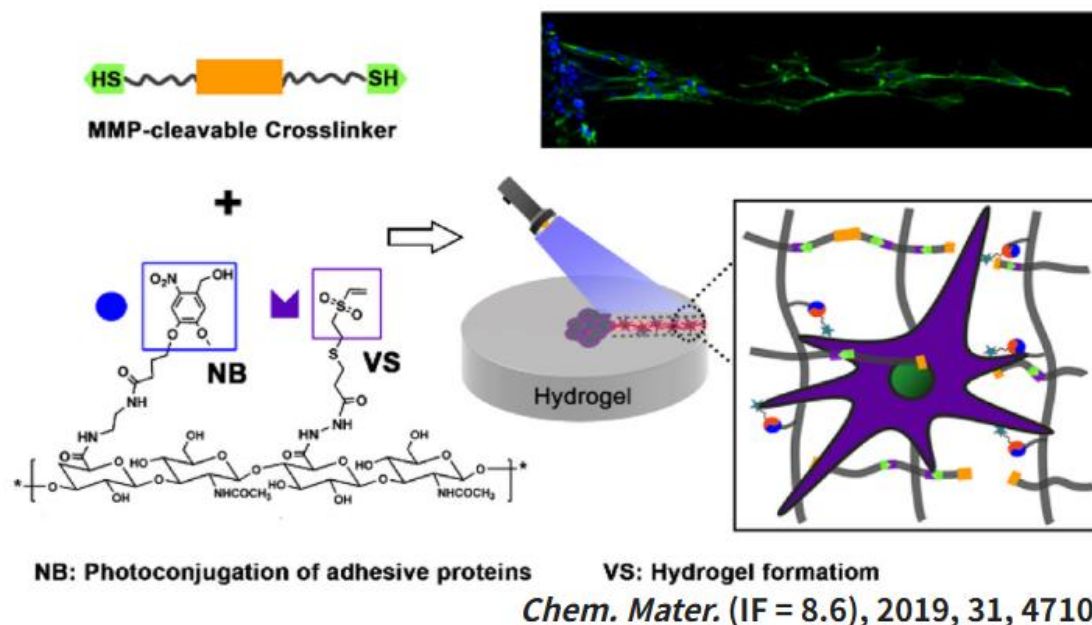


Fig. 9 Cell culture application.

Product Advantages

1. Batch-to-batch consistency: Standardized manufacturing ensures high repeatability and controllability in composition, physicochemical properties, and biological performance across batches.
2. Ultra-high purity: Unique purification process effectively removes residual NB monomers and process impurities, achieving an ultra-low impurity profile and ensuring safety in biomedical applications.
3. Reliable safety: HANB is the first commercially available product of its kind to have completed biocompatibility evaluation by a third-party institution, supporting safe and non-toxic use and accelerating lab-to-clinic translation.
4. High stability: In accordance with the Guidelines for Stability Studies of Non-Active Implantable Medical Devices, real-time and accelerated stability tests confirm that HANB remains stable under labeled storage conditions.

Specifications

Property	Value
CAS No.	N1518262

Property	Value
Molecular Weight	1180 kDa
Synonyms	o-Nitrobenzyl Alcohol-modified Sodium Hyaluronate HANB
Specifications & Purity	BioReagent, Molecular weight 1180 kDa; Degree of substitution 3.0%-4.0%; o-nitrobenzyl alcohol residue ≤1000 ppm
Stability And Storage	Store at -20 °C long term (24 months). Store in the dark.
Storage Conditions	Protected from light, Store at -20°C
Shipped In	Ice chest + Ice pads. This product requires cold chain shipping. Ground and other economy services are not available.

Contact & Global Offices

Whether you have a technical question, need help with a quotation, or want to inquire about an order, our regional teams are ready to assist. Please contact the office for your region; for general inquiries, the North American office is the corporate primary.

NORTH AMERICAN SALES, SUPPORT & GENERAL INQUIRIES

Aladdin Scientific Corporation

14078 Meridian Parkway, Riverside, CA 92518, USA

Phone: 1-833-552-7181

Sales: sales@aladdinsci.com

Customer Service: custserv@aladdinsci.com

EU SALES, LOGISTICS & LOCALIZED SUPPORT

Aladdin Biochem Deutschland GmbH

Westring 2, 33142 Büren, Germany

Phone: +02951 9383958

Support: support.eu@aladdinsci.com

Limitations & Disclaimer

- For Research Use Only (RUO). Not for use in human or animal diagnostics, therapeutics, or in vivo applications. Not for food, cosmetic, or household use.
- This product is not a CE-marked in vitro diagnostic device under IVDR (EU) 2017/746 and is not an FDA-cleared device under 21 CFR. Use is restricted to verified businesses, institutions, and qualified professionals for research and development purposes.
- Where any component is classified as hazardous under CLP (EC) 1272/2008 or OSHA HCS (29 CFR 1910.1200), the product Safety Data Sheet (SDS) takes precedence over this document for handling, storage, and disposal information.



ALADDIN SCIENTIFIC

Reagents · Kits · Standards

Product Instructions
For Research Use Only
www.aladdinsci.com

- Performance depends on sample type, sample condition, handling, and operator technique. Users are responsible for validating the product for their specific application.