

Kisser's Mounting Medium

K747449

Store at room temperature

Introduction:

The Kisser's Mounting Medium is an aqueous mounting solution prepared according to the classical Kisser method, which can be used for the mounting of various tissue sections and smears on slides to maintain samples in good condition. The main active ingredients contained in this product are gelatin and glycerin, ensuring a long-term storage of the Kisser's Mounting Medium and a stable preservation of samples at room temperature.

Usage method:

For Adherent cells:

- a. After staining and washing, remove and discard the wash solution.
- b. Add one drop of Kisser's Mounting Medium on the glass slide, cover with a coverslip with adherent cells allowing cells contact with the medium, and avoid the formation of air bubbles.

For tissue sections:

- a. After staining and washing, remove and discard the wash solution.
- b. Add one drop of Kisser's Mounting Medium onto the tissue section on a glass slide, cover with a coverslip and avoid the formation of air bubbles.

Precautions:

1. Kisser's Mounting Medium solidifies at low temperature. Thaw it completely at 37-50°C prior to use.
2. Kisser's Mounting Medium contains a small amount of phenol. Please be careful when handling the reagent and take effective measures to avoid inhalation and direct contact with skin or eyes.
3. This product is for R&D only. Not for drug, household, or other uses.
4. For your safety and health, please wear a lab coat and disposable gloves during the operation.